



Board Programming

Models: R1005 – R1050

Automatic 5L – Automatic 50L

To install replacement boards refer to the “RE-10 Series Board Replacement Guide”

The replacement boards we sent you are for a different model. They are all interchangeable with a very easy parameters update.

To begin, power on the repaired rotary evaporator.

Press and hold the silver knob until you see this appear on the screen:



Turn the knob until “68” is displayed:



Press the knob to accept.

Refer to the below table for your model. You will need to change values on your machine to reflect the parameters for your model.

<i>Parameter</i>	<i>Model RE-1005</i>	<i>Model RE-1020</i>	<i>Model RE-1050</i>
<i>"Cc"</i>	14	15.3	11.5
<i>"rP"</i>	12	12	24
<i>"Pr"</i>	5	5	5
<i>"2F"</i>	0	0	1
<i>"rA"</i>	4	4	4
<i>"m"</i>	120	120	120
<i>"P5"</i>	0000	0000	0000

Press the knob to move to the next parameter. On the last parameter "P5" Press the knob again to save the settings and close the parameter menu. You will be returned to the main menu.

Congratulations your Rotary Evaporator control board is programmed for your model.

If you have any questions or concerns, please give us a call at (734) 855-4890 or send me an email dan@usalabequipment.com