




RE-10 SERIES ROTARY EVAPORATOR MOTOR REPLACEMENT GUIDE

STEP 1: Remove the 4 screws on top of the motor housing.




STEP 2: Remove the 4 screws for the back plate.

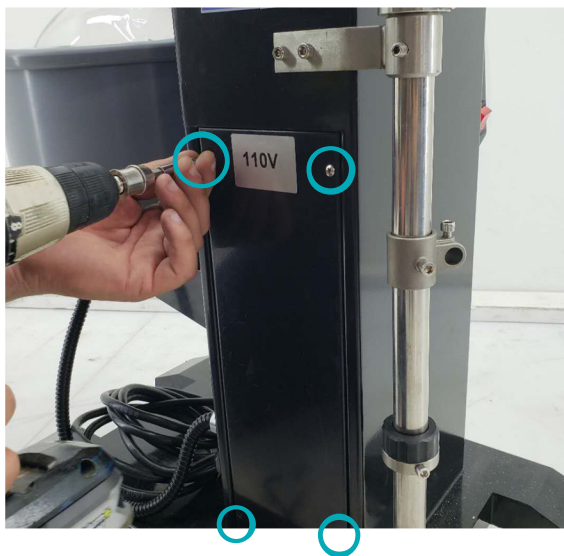


STEP 3: Cut the two zip-ties pictured below.

pg.2

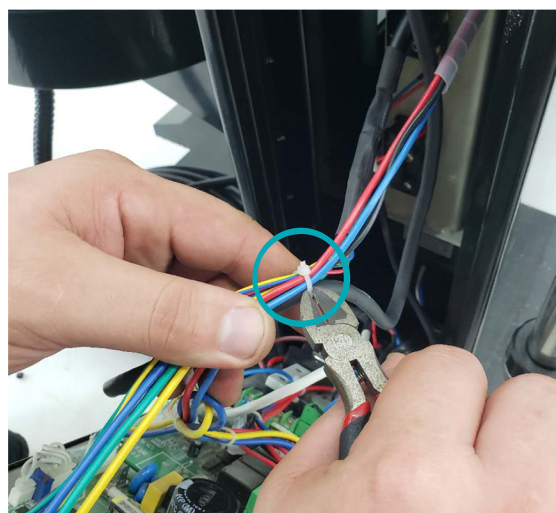

STEP 4: Disconnect the two connectors pictured below.

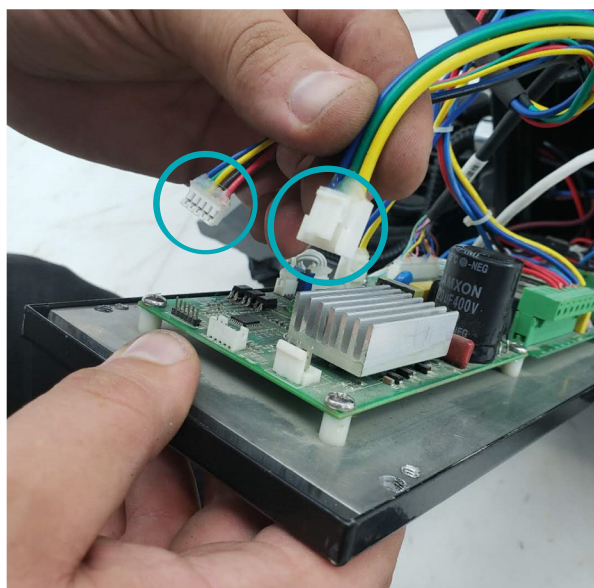

You can now remove the motor housing from the base. Once removed, you can re-install the motor and follow the steps in reverse.



If you have any questions or concerns, please sent us an email or give us a call at
734-855-4890 or sales@usalabequipment.com

THANK YOU!