



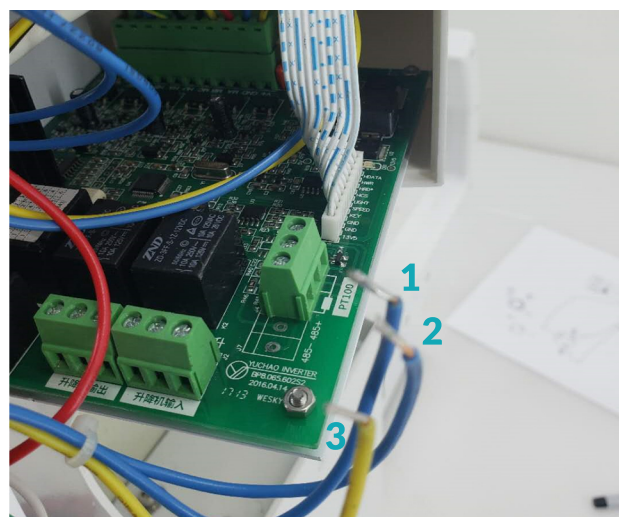


RE-200A ROTARY EVAPORATOR

PROBE CONNECTOR REPLACEMENT GUIDE

STEP 1: Once the back plate is removed, unscrew the 3 wires from the PCB connector.

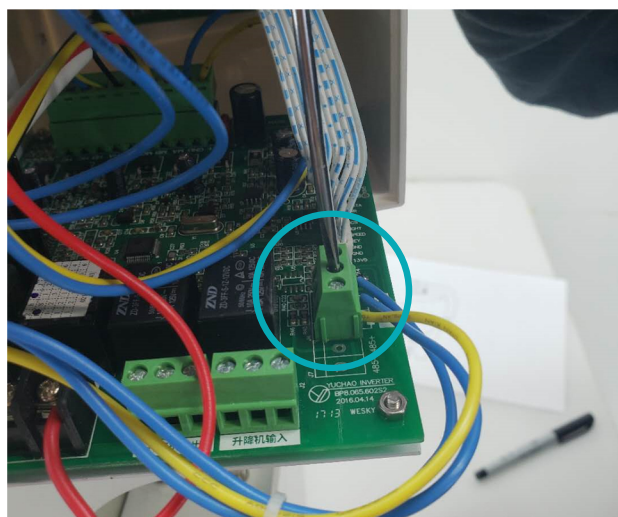


STEP 2: Unscrew the probe connector and remove it from the housing.

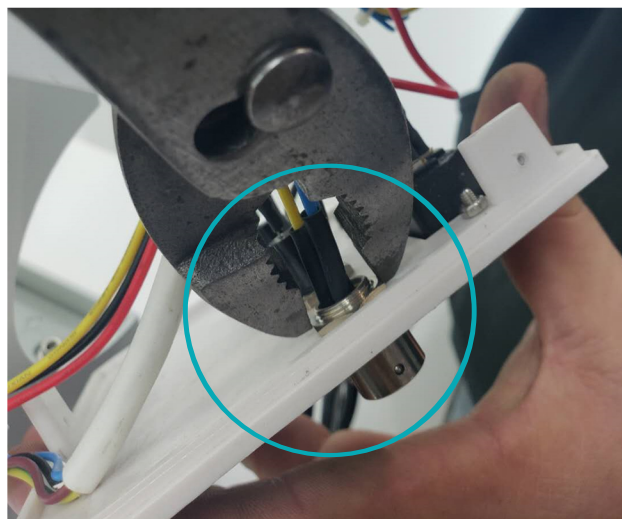


STEP 3: Insert the new probe connector into the housing and screw it on tightly.

pg.2

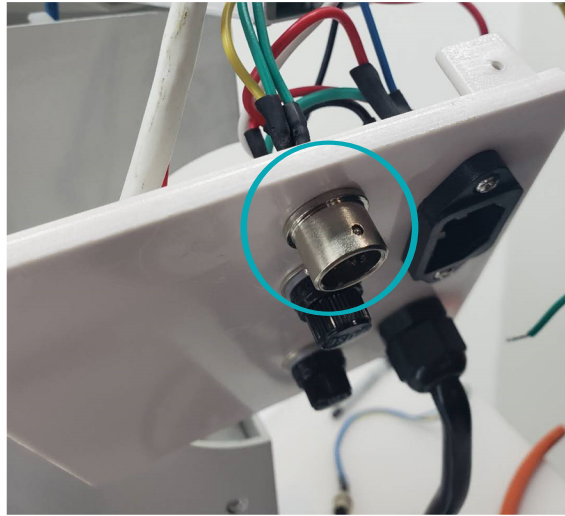


STEP 4: Screw all three wires into the PCB connector. (Wiring diagram below)

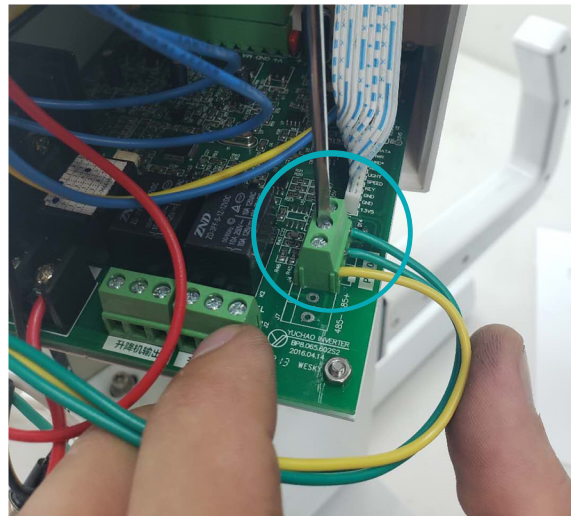

STEP 5: Re-connect all pieces and test the unit for proper operation.

WIRING DIAGRAM

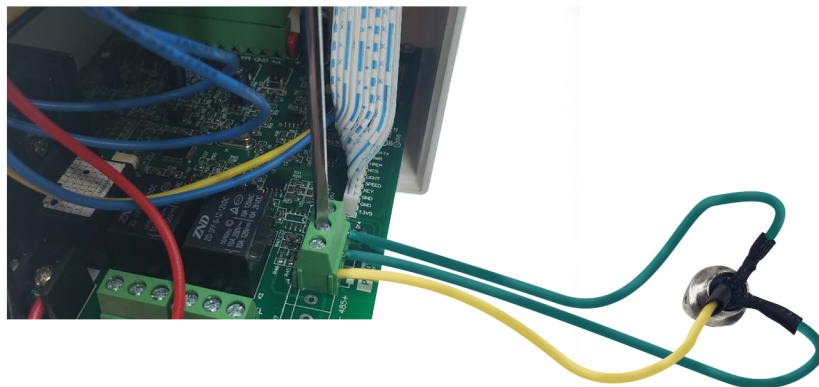

If you have any questions or concerns, please send us an email or give us a call at **734-855-4890** or sales@usalabequipment.com

THANK YOU!