




UC Series Pump Upgrade Instructions

(Includes HC-5/10)

[Step 1] Remove the old pump using our pump replacement guide.

- Resources tab on our website -

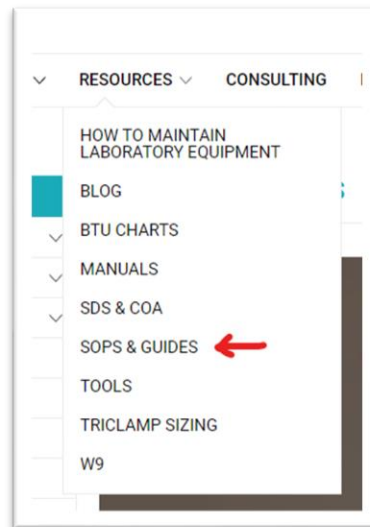

- Scroll to the file indicated below and open it -

• [USA Lab UC/DLSB-5/10 Series Chiller Pump Replacement Guide](#)


- Follow the guide steps 1-10 -

If you have any trouble locating our guide or need more information.


Please contact us.

tech@usalab.com | (734) 855-4890

[Step 2] Check the orientation of the new style pump vs. the old style.
(It may be necessary to remove the screws and rotate the housing 180° to match inlet and outlet path)

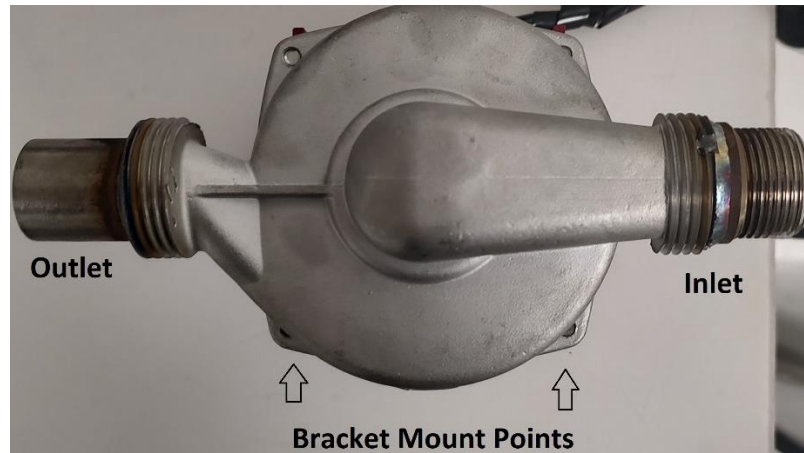

[Step 3] Attach pump to the old bracket by removing the two bottom screws. Add pump onto bracket and tighten screws.

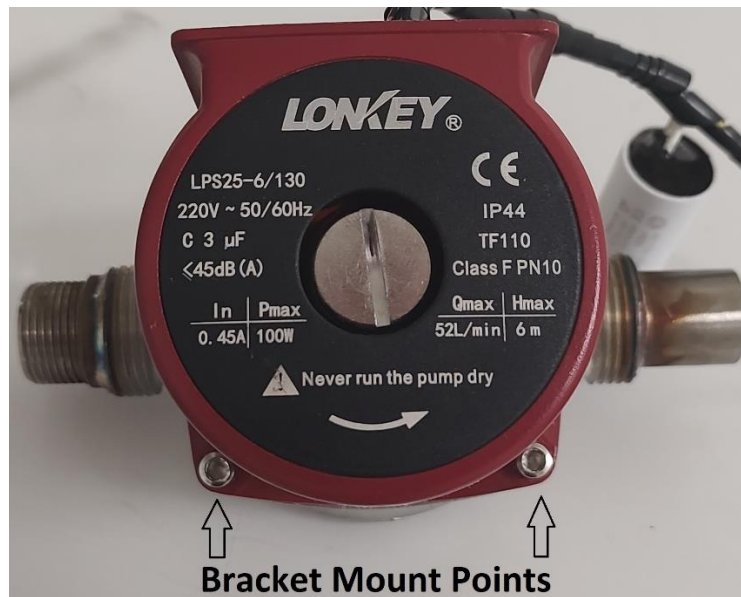

[Step 4] Re-wrap insulation onto the pump and reverse disassembly.

If you have any trouble or need more information.

Please contact us.

tech@usalab.com | (734) 855-4890