


# **OPERATION MANUAL Mobile Cooler**

MCF-1.09-25/MCF-2.82-25

**VERSION: 2017-4** 

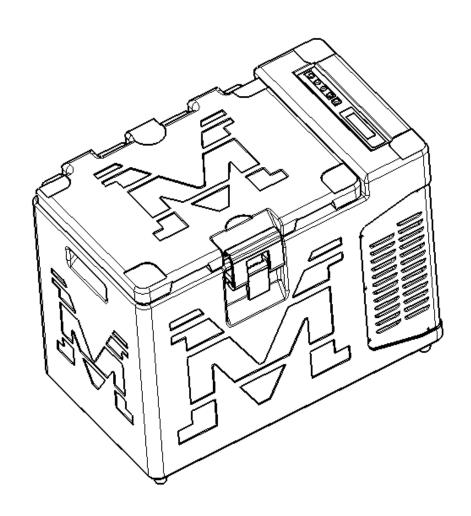



# **CONTENTS**

I. Introduction	2
1.1 Area of use	2
1.2 Consulting the manual	3
II. Warnings	3
2.1 General recommendations	3
III. Using the product	4
3.1 Using the product for the first time	4
3.2 Installation	4
3.3 Activation	6
3.4 Setup button	6
3.5 Loading instructions	7
3.6 Transport instructions	7
3.7 Instructions in case of prolonged disuse.	7
3.8 Cables	8
IV. Maintenance	8
4.1 Washing instructions	8
4.2 Defrosting instructions	8
4.3 Maintenance instructions	9
4.4 Solving problems	9
4.5 Replacing the protective fuse	11
V. Disposal	11
5.1 Packing	11
VI. Refrigerant gas	12
6.1 Product structure and composition	12
6.2 Main technical parameters	12
6.3 Circuit Diagram	13
VII. Packing list	13



## I. Introduction

## 1.1 Area of use

- > Thank you for your choosing our products ,For your safe, convenient use and reasonable maintenance of the products,Please carefully read the User's Manual before use and carefully store it for reference in the future.
- This series of medical vehicle low temperature storage cabinet have been designed and built for the transport of temperature sensitive foodstuffs, without interruption of the cold chain. They represent the solution for transport over long distances or for which precise temperature control is required for the duration of transport.
- ➤ Depending on the using condition, refrigerator can be divided into refrigerated transport (temperature control device is set to 0~+30 degrees Celsius) and frozen transport (temperature control device is set to -25~+30 degrees Celsius).
- ➤ According to different conditions, can be divided into refrigerated transport (thermostat setting range 0~+30 °C) and refrigeration transportation(thermostat setting range -25~+30 °C)
- ➤ The environment temperature of reference is between +10°C and +38°C. Outside of this range, the performance levels declared for the refrigerated containers may vary.

#### **Cautions on Safety**

- Must ensure that the storage refrigerator can not work in an environment of more than 38. If this limit is exceeded, the electronic part of the storage box will be completely destroyed.
- > Do not install the appliance in the damp site or place where may easily suffer from water, otherwise, it may cause electric leakage or electric shock since the insulating quality will be dropped.
- > Do not store flammable, explosive and dangerous products or volatile materials in the equipment, do not use flammable mist near the equipment, otherwise, it may possibly cause explosion or fire.
- > Do not store acid, alkali and corrosive materials in this equipment, otherwise, it may possibly damage the internal components or electric parts of the equipment.
- > Do not insert metal products such as iron nail or wire into any hole and clearance of this equipment or any outlet for internal air circulation, otherwise, it may possibly cause electric shock or physical injury since the above objects may contact with the moving parts.
- Please use the appliance in the safe area when it is used for toxic, harmful or radioactive materials since incorrect use may possibly cause harm to human health the environment.
- When there's flammable gas leakage, please turn off the gas valve, open the doors and windows for ventilation, please do not unplug or insert the power plug of the storage cabinet.
- Please cut off power supply before equipment repair or maintenance for avoiding electric shock or physical injury.
- > Ensure that the repair technician won't inhale medicine or suspended particles in or surrounding this equipment, otherwise, it may possibly cause harm to human health.
- Do not directly pour water on the appliance, otherwise, it may cause electric shock or short circuit.
- Please do not disassemble, repair or modify this appliance by yourself. Any above work which is not



- operated by the authorized technician may possibly cause fire or physical injury.
- The product is not sterilized, does not allow patients to contact, use or affect the product and product's storage;
- The product is only used for storage and it is not allowed to combined with other materials, organization and technology in the process of use.
- Products can not be used for measurement or analysis.
- The life of the product is not related to the life of storage goods.
- > Some models need to regularly add supplies, such as: printing paper, test tubes, etc.;
- > This product is not explosion-proof products
- Non frozen equipment, only use to maintain the temperature.

## 1.2 Consulting the manual

- > This manual has been arranged in such a way that the user can find all the information needed for quick and easy maintenance.
- Consequently it must be made available to maintenance staff and operators at all times.
- The instructions must be read carefully before using the product.

## II. Warnings

- > The product is built to the state of the art. All requirements for the safe and correct operation of the appliance have been satisfied.
- The manager of the user company is recommended to train staff in order to ensure that the container is destined exclusively to the use for which it has been designed and used correctly as indicated in these instructions.

#### 2.1 General recommendations

- > The use of the container must comply with the manufacturer's instructions. The function of the storage cabinet is that of guaranteeing the maintenance of temperature (chilled or frozen) required during transport.
- > Different functions or operating modes should be avoided.
- For repairs, always contact a service center authorized by the manufacturer and insist on the use of original spare parts.
- Never cover the ventilation grilles. Leave a free space of at least 20cm around the ventilation grilles.
- > Do not store the container into cold stores. The high humidity level in the air could damage the electronics of the container. The low temperature could solidify the oil contained in the refrigerated unit.
- > Do not store the container at a temperature lower than 0° C.
- > Periodically check for the condition of the container.
- > Use the container on flat surfaces only.
- > Switch off the cooling unit in the case of tipping or steep inclination. Re position the container on a flat surface and wait for at least 1 hour before switching the cooling unit on again.



- > Do not activate the cooling unit if it has been dropped or damaged.
- > Do not bring the container into contact with sharp or pointed surfaces or sources of direct heat (e.g.:naked flames).
- > Do not expose the container to splashing water, rain, weathering or aggressive and pollutant atmosphere (e.g:smoke, fumes).
- Production date see bar code, bar code first, second represent the year, Third, Fourth representatives batch number, fifth, sixth, seventh, eighth representatives of products units numbered, the last two on behalf of the production month.
- > Duration of use: Refer to the key components that affecting the use of the shortest life
- Never wash the container with steam jet or pressurized appliances.
- > Do not leave the container exposed to direct sunlight for a number of hours.
- > Do not put the storage box placed near the humidity or water leaching environment;
- > Do not install the container near cookers, heaters or other appliances which give off heat.
- > Please do not touch the internal of storage cabinet with a wet hand
- No corrosive, flammable, explosive gas, liquid or dust in the surrounding environment;
- > Ensure to keep the surroundings well ventilated.

The manufacturer declines all responsibility if the accident prevention standards in force are not observed by the user company.

## III. Using the product

## 3.1 Using the product for the first time

The container was cleaned before leaving the factory. However, we recommend washing the container, following the instructions given in paragraph 4.1 "Washing Instructions" before using it.

## 3.2 Installation

- > This series refrigerator can be powered with 12V DC power or with 110V AC 50/60Hz. power.
- For the DC use, This product can be applied to a battery-powered motor vehicles for the use of 12V.
- For the AC use, the refrigerator are set for standard operation with 240V AC; for the operation with 100V AC it is necessary to operate a selector located on the power supply(part of the product can only be applied 110V).

If the local grid voltage is too high or too low, the compressor does not work and the electronics of the equipment may be damaged.

The storage cabinet is protected by the following fuses:

- > 15 A fuse, located near the 12V DC socket for the refrigerator with integrated cooling unit
- > 12A fuse, located on the internal power supply 12V DC cigarette lighter

#### Alternate current connection 110V AC 50/60Hz:

- > check that the plug on the power cable is suitable for the electricity socket;
- ensure that the socket has an efficient earth contract and adequate capacity. The electrical safety is only



guaranteed when it is correctly connected to an efficient earth system; systems which do not comply with the standards in force could cause injury or damage;

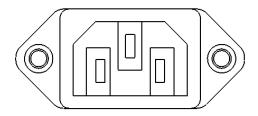


Figure 1:110 v ac 50/60 HZ connection socket and fuse

#### 12V DC connection:

- use original cables supplied by the manufacturer only;
- different connections must be assessed and made by qualified personnel;
- in case of installation of the refrigerator on the same vehicle, We recommend to contact manufacturers after-sales department, in order to evaluate the correct current inputs and the sections of the power cables, to guarantee the correct operation of the system.

#### To prevent drops in voltage and losses of power:

- The cable must be as short as possible and must not be sectioned;
- Avoid using additional switches, plugs or junction boxes;
- Do not connect other electric appliances on the cable of the cooling unit;
- > The connection of the cooling unit to the battery must be direct and dedicated;
- Do not use portable generators, they could cause damages due to voltage peaks and frequency variations;
- Do not use battery chargers to power the cooling unit;
- Maintain a free surface around the cooling unit (at least 20 cm), to ensure an adequate ventilation and allow a greater cooling efficiency and a reduced power consumption.

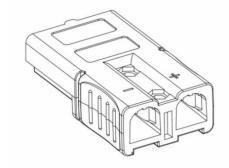


Figure 2: Refrigerator 12V DC connection socket



## 3.3 Activation

#### 110V 50/60Hz

#### 50/60Hz 110V AC connection

- push the power cable socket into the appliance socket (Drawing 1);
- > push the mains plug into the electric socket 110V AC 50/60Hz;
- The main switch (O/I) is pushed to the position of I (O/I), storage cabinet work (see Drawing 3);

#### 12V DC link

- push the power cable socket into the appliance socket (Drawing 2);
- connect the cooling unit to the 12 V DC power supply;
- Push the main switch (O/I) to the position I, the storage cabinet start working (see Figure 3).
- > according to the storage cabinet model, The temperature regulation device is programmed in the factory at a set-point value.

## 3.4 Setup button

- > There are six separate buttons 🗸 🔯 🔨 🗸 🐧 on the control panel for control of the thermostat;
- → + A Keyboard lock; A + Defrost control; Ø + A + Reset; A + A C setting; A + A 22 C setting.

#### **Description** Power Switch

- ➤ Main Power open: Press to control the display of the controller;
- Main Power open: Press 0 ≥6s can turn on or off refrigeration system related components;
- Factory default temperature display and refrigeration system related components are open state.

#### Setting button

- In the normal mode, press ⇒ ≥1s, will be enter the temperature modification mode, you can change the setting temperature;
- In the temperature modification mode, press to save the settings and return to the normal display state, otherwise automatically give up the changes after 6s and return to the normal display state;

## △Up

➤ When setting temperature, add set temperature value by press :

### **∨** Down

➤ When setting the temperature, press , reduce the set temperature value.

#### Mute button

Press \( \) to stop buzzer.





**Drawing 4: Temperature controller panel** 

#### Alarm code

Alarm code	Alarm description	Alarm code	Alarm description
АН	Cabinet temperature too high alarm	AS1S	Control probe 1 short circuit
AL	Cabinet temperature too low alarm	AS1B	Control probe 1 short circuit
AS2S	Control probe 2 short circuit	AS2B	Control probe 2 short circuit
AUF	Power failure alarm		

## 3.5 Loading instructions

- Open the storage cabinet using the plastic handles.
- The container is suitable for contact with items; this means that loose or unpacked food can be stored inside of it.
- > The purpose of storage cabinet is not to heat or cool the products, but to guarantee that they maintain their temperature. Therefore it is vital to load the goods inside the container at the temperature required.
- ➤ It is advisable to arrange the products inside without cardboard packaging as this delays the penetration of the cold.
- To avoid the dispersion of heat, keep the container open as briefly as possible.
- Close the container using the special locking handles immediately after storage the items.

## 3.6 Transport instructions

- Before moving the container, make sure that the lid or the door is closed.
- The loaded storage cabinet may reach a considerable weight; Therefore must always be careful to move. If possible, you need to use special car. Ascension, transportation, storage cabinet, should be used to save the sag of both side of the type of handle
- Use the appropriate handles to move the container provided with wheels or trolley.
- When moving the container, pay the utmost attention to avoiding collision with objects or people.
- To avoid storage cabinet door open in the ordinary state, it may causing the temperature rise rapidly.

## 3.7 Instructions in case of prolonged disuse.

- If the container is not used for a prolonged period of time, carry out the following operations:
- remove all the products from the container;
- > switch off the cooling unit by pressing 0 on the main switch O/I (Drawing 3). Pull out the plug connected to the electricity main (or battery) and the plug connected to the appliance. Store the cable in a safe place and protect it against damp;



- clean the container as indicated in paragraph 4.1 "Washing instructions";
- leave the lid (door) open for a few hours to prevent the formation of unpleasant smells.

#### 3.8 Cables

- We recommend the use of original cables only. Different connections must be assessed and made by qualified personnel.
- > To prevent drops in voltage and losses of power, the cable must be as short as possible and must not be sectioned; therefore, it is necessary to avoid using additional switches, plugs or junction boxes.
- > The section of the cable must be chosen in accordance with its length.
- ➤ In particular:

### For connection to the 12V DC power supply:

2P Cable 1.8 metres long, section 2.5 mm², for storage cabinet .

#### For connection to the 110V AC 50/60Hz power supply:

- > 3P Cable 2 meters long, section 0.75 mm<sup>2</sup>.
- In case of installation of multiple storage cabinet on the same vehicle, we recommend to refer to the after-sale department of the manufacturer, in order to evaluate the correct current inputs and the sections of the power cables, to guarantee the correct operation of the system.

## IV. Maintenance

## 4.1 Washing instructions

- Regularly clean the container inside and out, observing the following instructions:
- Before cleaning the container, switch off the cooling unit by pressing 0 on the main switch O/I (Drawing 3). Pull out the plug connected to the electricity main (or battery) and the plug connected to the appliance. Store the cable in a safe place and protect it against damp.
- ➤ Clean the container using a cloth and warm water; for the purposes of hygiene, always dry the container using disposable cloths or paper (never reusable cloths).
- > Do not use metal or synthetic scouring pads, only soft brushes with natural or plastic bristles.
- > Do not use abrasive powder, ammonia, acid or solvent.
- It is possible to use soap solutions.
- Do not use steam jet or pressurized appliances.
- Ensure that the water does not penetrate the power and regulation controls, vents or appliance socket.

## 4.2 Defrosting instructions

- If the external temperature and humidity are high and the lid (door) is opened frequently, a layer of frost will gradually develop on the surface of the evaporator. This layer acts as insulation and, if it exceeds a thickness of 3 mm, may reduce cooling efficiency. This makes regular defrosting necessary, as follows:
- remove all the products from the container;
- > switch off the cooling unit by pressing 0 on the main switch O/I (Drawing 3). Pull out the plug connected



- to the electricity main (or battery) and the plug connected to the appliance. Store the cable in a safe place and protect it against damp;
- > leave the lid (or door) of the container open, until the layer of frost has completely melted. Do not try to speed up the defrosting process using heaters or remove frost using knives or other sharp objects;
- After the defrosting, dry the container using a disposable cloth or paper. The container is now ready to be used.

## 4.3 Maintenance instructions

Before carrying out any maintenance operation, switch off the cooling unit by pressing 0 on the main switch O/I (Drawing 4 - A). Pull out the plug connected to the electricity main (or battery) and the plug connected to the appliance. Store the cable in a safe place and protect it against damp.

#### We recommend:

- > maintenance of the cooling unit, carried out by qualified technicians at least once a year;
- use of original spare parts at all times.
- > Do not allow children to use medical freezer to playing games .
- Not to put heavy objects on the refrigerator door or medical cap, to avoid the door or the top compression deformation.
- Contact your local dealer or the after-sale department of the manufacturer for assistance and supply of original spare parts.

#### **ORDINARY MAINTENANCE**

PART TO CHECK	FREQUENCY	TYPE OF CHECK
Power cable and main plug		Check that they are neither damaged nor too
Fower cable and main plug	6 months	old.Replace them if necessary.
		Check the condition.
Seal	6 months	Replace it if broken or worn.
		frost is more than 3 mm thick, defrost (following
frost on the surface of the evaporator	One week	the instructions given in paragraph 4.2
	One week	"Defrosting instructions")

## 4.4 Solving problems

The table below lists the major faults that can be found on the refrigeration unit of the storage cabinet, with indication of the possible causes and interventions for recovery.

If you can't judge fault reason, or difficult to exclude failure, please according to the warranty certificate (random annex) instructions and we stay in a warranty certificate after-sales contact, inform your name, address, telephone number, for medical refrigerator types, purchase time, fault phenomena, the professional engineer will wholeheartedly at your service.

FAULT	POSSIBLE CAUSE	TYPE OF INTERVENTION
The cooling unit does not operate with a 12V DC power supply	a) The refrigerator is not connected to the 12V DC power supply	Connect the cooling unit to the 12V DC power supply. Check the 12V DC power cable and replace it if necessary
	b) The 12V DC line fuse	b) Install a new fuse on the 12V DC line. (the



	has blown	fuse is
	TIAS DIOWIT	located near the 12V DC socket (Paragraph
		4.6).
		Investigate the cause of the failure (likely
		short circuit or momentary voltage overload).
	c) The battery is dead	c) Test the battery and charge it or replace it.
	d) The main switch 0/I	d) Check the main switch 0/I (Drawing 3)
	(Drawing 4 - A) is faulty	and replace it if necessary.
	e) The electric wiring is	e) Restore the wiring as shown in the wiring
The cooling unit does not	disconnected	diagram.
The cooling unit does not operate with AC power supply	a) The cooling unit is not connected to the 110V AC	Connect the cooling unit 110V AC 50/60Hz power supply.
operate with AC power supply	50/60Hz power supply.	Check the 110 V AC
	oo, oor 12 power suppry.	50/60Hz power cable and replace it if
		necessary.
	b) The fuse of the 110V AC	b) Install a new fuse on the power line and
	50/60Hz line is faulty	check the failure reason (likely short-circuit
		or momentary voltage overload).
	c) The main switch 0/I	c) Check the main switch 0/I (Drawing 3)
	(Drawing 3) is faulty.	and replace it if necessary.
	d) The electric wiring is	d) Restore the wiring as shown in the wiring
The coolings with door and	disconnected	diagram
The cooling unit does not maintain the set temperature	a) The ventilation is not sufficient for the cooling	a) Check that the ventilation grilles are free.
	b) The room temperature is	b) Let air circulate in the area where the
	too high	cooling unit is located.
	c) The evaporator is	c) Defrost as indicated in paragraph 4.2.
	covered in frost	
	d) The products are stored	d) Cool the products before putting them in the container
	at an excessively high temperature	the container
	e) The lid (or door) does	e) Check lid (or door)closing and possibly
	not close correctly	replace the seal.
	f) The temperature	f) Check the setting of the temperature
	regulator is not set	regulator and change the set-point value if
	correctly	necessary.
	g) The cooling system has failed	g) Contact the service center.
The temperature regulator is	a) The environment is too	a) Cover the temperature regulator
not turned on	bright	
	b) The main power	b) Insert the plugs into the right sockets.
	cable/battery is not	
	connected	
	c) The main power cable	c) Check the power cable to the grid/battery
	/battery is damaged d) The temperature	<ul><li>and replace it if necessary.</li><li>d) Check the temperature regulator and</li></ul>
	regulator electronics have	replace it if necessary.
	failed	replace it il fleecosary.
The temperature regulator	a) Faulty electrical contact	a) Check the correct supply of the
flashes continuously	, , , , , , , , , , , , , , , , , , , ,	temperature regulator.
	b) The temperature	b)replace
	regulator is faulty	
The temperature regulator	a) The temperature probe	a) Check the temperature probe and replace
displays numbers outside of	is faulty	it if necessary
the standard numbering	b) The connection of the	b) Check the connection of the probe with
	probe with the temperature	the regulator and restore it if needed.
The temperature regulator is	regulator is not correct  a) The pickup voltage is	a) The battery is dead: recharge or replace it
turned on but the compressor	lower than 11V	if necessary
tarried on put the compressor	I TO WOLL II II II II II	ii nooosai y



1 1 1 1 10 1 00		1) 01 1 4 1 1 4
does not start with 12V DC	b) The cable sections are	b) Check the power cables and replace them
power supply	not correct (inadequate	if necessary
	extensions)	
	c) The temperature	c) Check the setting of the temperature
	regulator is not set	regulator and change the set-point value if
	correctly	necessary (see paragraph 3.4 "Entry into
	-	service").
The temperature regulator is	a) The pickup voltage is in	a) The battery is almost drained: recharge or
turned on but the compressor	the range 11V to 11.5V	replace it if necessary.
does not always start with	b) The cable sections are	b) Check the power cables and replace them
12/24V DC power supply	not correct (inadequate	if necessary
	extensions)	
	c) The room temperature is	c) Let air circulate in the area where the
	too high	cooling unit is located.
The refrigeration unit drops in	a) Partial leak of gas	a)A) check the pressure and temperature in
temperature only by a few	a, raman ram or gas	refrigeration loop (contact after-sales
degrees, but works		personnel and local maintenance network to
degrees, but works		determine the location of the leak and
		supplementary refrigerant)
	b)Insufficient supply	b) Ensure that the power supply is suitable.
	voltage	b) Ensure that the power supply is suitable.
The refrigeration units noisy	a) Noisy fans	a) Check the operating status of the fans
and has strong vibration	a) Noisy lans	and replace them if necessary.
and has strong vibration	b) Motor unit not proporty	
	b) Motor unit not properly	b) Check the correct mounting of the motor
T1	fixed	unit and restore it if needed.
The refrigeration unit falls		Position the container on a flat surface.
over and overturns		Open the cover of the refrigeration unit,
		inspect the status of the components, check
		the pipes and electrical
		connections. If you are having no special
		problems, wait at least 1 hour before
		restarting the refrigeration unit.

## 4.5 Replacing the protective fuse

#### To replace the fuse on the 12V DC line, proceed as follows:

- > pull out the plug connected to the 12V DC power line and the plug connected to the appliance;
- open the fuse box using a tool (e.g.: the tip of a screw driver);
- replace the fuse and insert it into the fuse holder;
- plug in the plug connected to the appliance and the socket to the 12V DC power line.

# V. Disposal

## 5.1 Packing

The packing material (cardboard, polyethylene film) is 100% recyclable.

The user is responsible for disposal in observance of the local regulations.



# VI. Refrigerant gas

- > The cooling gas HFC 134a used in the cooling circuit of the refrigerator is not noxious-to the ozone
- > The cooling circuit is hermetic, without the possible leakage of coolant under normal conditions of operation and use.
- > The compressor used is hermetic, designed for applications in motion.

## 6.1 Product structure and composition


This product is mainly composed of rotational body, a door body, a refrigeration system, control system and accessories.

## 6.2 Main technical parameters

Model	Climate type	Temperature in Cabinet (℃)	Effective Capacity ( L )	Rated voltage $(V\sim)$	Refrigerant	External size (W*D*H)
MCF-1.09-25	SN/N	<b>-</b> 10∼-25	30	DC12/ AC110V	R134a	615*370*471.5mm
MCF-2.82-25	SN/N	<b>-10∼-25</b>	80	DC12/ AC110V	R134a	937*565*580.5mm



## **6.3 Circuit Diagram**

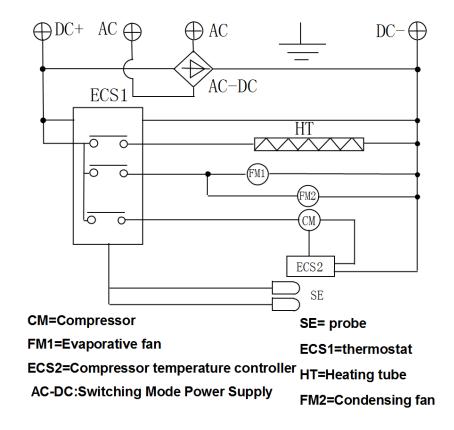

# VII. Packing list

Model	Manual	12V DC power line	110V AC power line	Plastic frames
MCF-1.09-25	1	1	1	1
MCF-2.82-25	1	1	1	1



(	Certificate of Quality	
Inspector:		