


Ultrathin Magnetic Stirrer Flat spin

Safety Warning

- The product should be operated only by trained staff only.
- Please do not operate this instrument in any manner not mentioned in this manual.
- If the instrument gives any strange noise or other abnormal phenomena while in operation, please immediately shut down the instrument and contact our service center.
- Keep away from the effects of high magnetic fields.
- Ensure that there is no possibility for any liquid spillage.

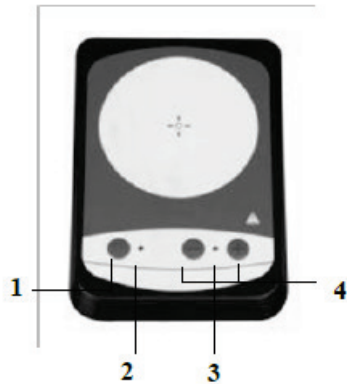

Operating Instructions

Switch ON/OFF:

- 1) **Switch ON:** Long press button 1 (power button) (for more than 1sec.) to switch **ON** the instrument, function indicator 2 lights up.
- 2) **Switch OFF:** Long press the button 1 (power button) (for more than 1sec.) to switch **OFF** the instrument, both indicators turn off.

Turn ON Forward/Reverse Mode:

Press button 1 (power button) to switch ON or OFF the reverse mode and invert the stirring direction at regular intervals.

- 1) **On state:** The stirring direction will be inverted every 15 seconds. The function indicator 2 blinks when the reverse mode is active.
- 2) **Off state:** The stirring direction will stay the same, function indicator 2 keeps lighting up.

Speed setting:

- 1) Press button 4 (speed adjustment button) to reduce or increase the speed, the speed setting is at intervals of 10rpm.
- 2) Long press the button 4 (speed adjustment button) to change the set speed continuously, the speed indicator 3 flashes frequently during the operation.
- 3) When the speed is set to a max. of 1500rpm or min. speed of 15rpm, the display indicator 3 will stop flashing.
- 4) The display indicator 3 will stop flashing and keep bright when the actual speed reaches to set speed value.

Common faults:

When the stirring bar supplied with this instrument does not work or any key malfunctions, Please connect with direct dealer.

12301837



Please scan the QR code to get more instructions